

Cas13 [LwCas13a (LwC2c2)] Kit

Catalog # CS13A-H834

Product Description

LwCas13a enzyme reaction kit consisting of enzyme, buffer, and optional reporter and/or guide RNA.

Components

	Component Name	20 µg	100 µg
CS13A-E321H	LwCas13a, Active	20 µg	100 µg
CS13-09B	Cas13 Reaction Buffer II, 10X	1ml	2 x 1ml
CS13-57	Cas13 ssRNA Reporter (optional)	10 nmol	10 nmol
CS13-G58	Cas13 Guide RNA (optional, custom)	1 vial	1 vial

Formulations

Enzyme	50 mM Tris-HCL pH7.5, 600 mM NaCl, 2 mM DTT, 10% glycerol
Buffer	200mM HEPES, pH 7.0, 500mM KCl, 50mM MgCl ₂ , 1mg/mL Prionex® (0.1% w/v)
Reporter	Lyophilized powder
Guide RNA	Nuclease free water

Storage and Stability

Store kit at -70°C. To avoid repeated handling and multiple freeze/thaw cycles aliquot components into smaller quantities. Protect reporter component from light.

Always include RNase inhibitor at manufacturer recommended concentrations in Cas enzyme reactions.

Scientific Background

Cas13a is an RNA-guided RNA endonuclease that belongs to the class 2 type VI CRISPR-Cas system (1). Cas13a possesses both RNA recognition and cleavage activities and cleaves RNA through conserved basic residues within its two HEPN domains (2). Activated Cas13a also induces collateral cleavage of nearby non-targeted RNAs in a nonspecific manner (3). This property of the Cas13a enzyme has been leveraged for the development of sensitive and robust nucleic acid detection tools with applications in both basic research and therapeutics development (4).

References

1. East-Seletsky A, et al. 2016. Two distinct RNase activities of CRISPR-C2c2 enable guide-RNA processing and RNA detection. *Nature*. 538:270–3.
2. Abudayyeh O, et al. 2016. C2c2 is a single-component programmable RNA-guided RNA-targeting CRISPR effector. *Science*. 353:5573.
3. Zhang J, You Y. 2020. CRISPR-Cas13a system: A novel approach to precision oncology. *Cancer Biol Med*. 17(1):6–8.
4. Patchesung M, et al. 2020. Clinical validation of a Cas13-based assay for the detection of SARS-CoV-2 RNA. *Nat Biomed Eng*. 4(12):1140-1149.

Reaction Protocol

Other Materials Required

- Filter pipette tips
- Nuclease free water
- Nuclease free microcentrifuge tubes
- Half-area solid black 96-well plate
- Microplate sealing tape
- Fluorescent microplate reader
- ssRNA substrate

Step 1: Thaw the active enzyme on ice. Prepare 1X Reaction Buffer with nuclease free water. Reconstitute reporter in nuclease free water. Equilibrate the buffer, reporter, guide RNA, and ssRNA substrate to ambient temperature.

Step 2: Prepare the following working solutions with 1X Reaction Buffer:

- 4X final concentration of Active Cas13 enzyme
- 4X final concentration of guide RNA
- 2X final concentration of substrate/reporter mix

Step 3: In a half-area solid black 96-well plate, add the following components and pre-incubate at room temperature for 10 minutes;

Component 1. 10 µL of 4X Active Cas13

Component 2. 10 µL of 4X guide RNA

Note: A blank control can be set up as outlined in step 3 by replacing the enzyme working solution with an equal volume of the reaction buffer.

Step 4: To each assay well, add 20 µL of the 2X substrate/reporter mix. Shake the plate for 1 minute on a tabletop orbital shaker. Seal the assay wells with microplate sealing tape and incubate at 37°C for 15-30 minutes.

Step 5: Equilibrate the plate to ambient temperature and then remove the microplate sealing tape. Read fluorescence on a microplate reader.

To place your order, please contact us by phone 1-778-326-0223 or 1-888-606-3424 (Toll free) or by email: orders@signalchemdx.com or info@signalchemdx.com - www.signalchemdx.com

FOR IN VITRO RESEARCH PURPOSES ONLY. NOT INTENDED FOR USE IN HUMANS OR ANIMALS.

SAFETY DATA SHEET

Article 1 - Product Identification

Product Name: Cas13 [LwCas13a (LwC2c2)] Kit

Catalog # CS13A-H834

This product is sold only for research use by qualified laboratory personnel, and is not to be used as a drug, medical device, food additive, cosmetic, nor household chemical. It is not to be used in diagnostic, therapeutic, consumer, agricultural, nor pesticidal applications.

Supplier of Datasheet: SignalChem Diagnostics Inc.
 Street Address: 190-13160 Vanier Place
 City, Prov. Postal Code: Richmond, BC, V6V 2J2
 Country: Canada
 Emergency Phone: 1-888-606-3424 (Toll free)
 1-778-326-0223 (local)

Article 2 - Hazard Identification

- **Emergency Overview:** The product contains no substances which, at their given concentrations, are considered to be hazardous to health.
- **WHMIS Classification:** Not WHMIS controlled.
- **GHS classification:** Not classified.
- **Hazard Pictograms:** None.
- **Signal words:** None.
- **Hazard statements:** None.
- **Precautionary statements:** None.
- **Other hazards:** None known.

Article 3 – Composition/Information on Ingredients

Description: This product consists of the components listed below

Component: LwCas13a, Active

Chemical Characterization: Mixture.

Common name	Chemical name	CAS-No.	Concentration
Glycerol	Glycerol	56-81-5	10%
NaCl	Sodium chloride	7647-14-5	3.5%
Tris	Tris(hydroxymethyl)aminomethane	77-86-1	0.61%
DTT	1,4-Dithio-DL-threitol	3483-12-3	0.0309%
Protein	N/A	N/A	≤0.01%

Component: Cas13 Reaction Buffer II, 10X

Chemical Characterization: Mixture.

Common name	Chemical name	CAS-No.	Concentration
HEPES	N-(2-Hydroxyethyl)piperazine-N'-(2-ethanesulfonic acid)	7365-45-9	4.8%
KCl	Potassium Chloride	7447-40-7	3.7%
MgCl ₂	Magnesium Chloride Hexahydrate	7791-18-6	1%
Prionex®	Prionex® Reagent	N/A	0.1%

Component: Cas13 ssRNA Reporter (optional)

Not a hazardous substance or mixture

Component: Guide RNA (optional custom)

Not a hazardous substance or mixture

To place your order, please contact us by phone 1-(604)-232-4600, fax 1-604-232-4601 or by email: orders@signalchem.com or IVD@signalchem.com - www.signalchem.com

FOR IN VITRO RESEARCH PURPOSES ONLY. NOT INTENDED FOR USE IN HUMAN OR ANIMALS.

SAFETY DATA SHEET

Article 4 – First-aid Measures

- **General information:** Consult a physician by providing the SDS.
- **After inhalation:** Breathe in fresh air. If casualty cannot breathe, give artificial respiration and consult a physician.
- **After skin contact:** Immediately wash with soap and plenty of water and rinse thoroughly. Consult a physician.
- **After eye contact:** Rinse opened eyes with plenty of water for at least 15 minutes. Remove contact lenses, if present and easy to do so. Consult a physician.
- **After swallowing:** Not expected to present a significant ingestion hazard under anticipated conditions of normal use. If you feel unwell, seek medical advice.

Article 5 - Fire-fighting Measures

- **Suitable extinguishing media:** Use water spray, extinguishing powder, carbon dioxide, or other appropriate measure that is suitable to the environment.
- **Specific hazards arising from the substance or mixture:** None known.
- **Special protective equipment and precautions for fire-fighters:** Self-contained breathing apparatus if necessary.

Article 6 – Accidental Release Measures

- **Personal precautions, protective equipment and emergency procedures:** Apply standard laboratory practices and personal protective equipment. Avoid breathing vapors, mist, or gas. Ensure adequate ventilation.
- **Environmental precautions:** Do not allow to enter drains.
- **Methods and materials for containment and cleaning up:** Absorb on sand or vermiculite and place in closed containers for disposal.

Article 7 - Handling and Storage

- **Precautions for safe handling:** Wear chemical safety goggles and compatible chemical-resistant gloves. Avoid inhalation, contact with eyes, skin or clothing.
- **Conditions for safe storage:** Store according to product label instructions. Keep container upright and tightly closed.

Article 8 - Exposure Controls/Personal Protection

- **Components with limit monitoring values at workplace:**
NA
- **Appropriate engineering controls:**
Apply adequate ventilation including mechanical exhaust or laboratory fume hood. Follow standard laboratory practices.
- **Individual protection measures:**
Respiratory protection:
Use appropriate respirator if there is inadequate ventilation by following the government standards.
Hand protection:
Wear gloves and use proper glove removal technique to avoid skin contact. Discard gloves after use by following the applicable laboratory regulations. Wash and dry hands.
Eye/face protection:
Safety goggles with side-shields approved under appropriate government standards.
Skin/body protection:
Use appropriate clothing, footwear and any additional protection measures to protect from splashing or contamination.

Article 9 – Physical and Chemical Properties

Component: LwCas13a, Active

Appearance: Colorless fluid.	Danger of explosion: Product does not present an explosion hazard.
Odour/Odour Threshold: Not determined.	Explosion limits: Not available.
pH: Not available.	Decomposition temperature: Not available.
Melting point/freezing point: Not determined.	Vapor pressure at 20 °C: Not available.
Boiling point/Boiling range: Not determined.	Density: Not determined.
Flash point: Not determined.	Relative density: Not determined.
Flammability (solid, gaseous): Not determined.	Vapor density: Not determined.
Ignition temperature: Not determined.	Evaporation rate: Not determined.
Auto-igniting: Product is not self-igniting.	Solubility in / Miscibility with Water: Fully miscible.

To place your order, please contact us by phone 1-(604)-232-4600, fax 1-604-232-4601 or by email: orders@signalchem.com or IVD@signalchem.com - www.signalchem.com

FOR IN VITRO RESEARCH PURPOSES ONLY. NOT INTENDED FOR USE IN HUMAN OR ANIMALS.

SAFETY DATA SHEET

Component: Cas13 Reaction Buffer II, 10X

Appearance: Colorless fluid.	Danger of explosion: Not determined.
Odour/Odour Threshold: Not determined.	Explosion limits: Not available.
pH: Not available.	Decomposition temperature: Not available.
Melting point/freezing point: Not determined.	Vapor pressure at 20 °C: Not available.
Boiling point/Boiling range: Not determined.	Density: Not determined.
Flash point: Not determined.	Relative density: Not determined.
Flammability (solid, gaseous): Not determined.	Vapor density: Not determined.
Ignition temperature: Not determined.	Evaporation rate: Not determined.
Auto-igniting: Not determined.	Solubility in / Miscibility with Water: Fully miscible.

Component: Cas13 ssRNA Reporter (optional)

Appearance: Translucent or opaque flakes, film or powder like substance	Danger of explosion: Not determined.
Odour/Odour Threshold: Not determined.	Explosion limits: Not available.
pH: Not available.	Decomposition temperature: Not available.
Melting point/freezing point: Not determined.	Vapor pressure at 20 °C: Not available.
Boiling point/Boiling range: Not determined.	Density: Not determined.
Flash point: Not determined.	Relative density: Not determined.
Flammability (solid, gaseous): Not determined.	Vapor density: Not determined.
Ignition temperature: Not determined.	Evaporation rate: Not determined.
Auto-igniting: Not determined.	Solubility in / Miscibility with Water: Fully soluble.

Article 10 - Stability and Reactivity

- **Reactivity:** Stable under recommended transport and storage conditions.
- **Chemical stability:** Stable under recommended transport and storage conditions.
- **Possible hazardous reactions:** No dangerous reactions known.
- **Conditions to avoid:** Heat and moisture.
- **Incompatible materials:** Not determined.
- **Hazardous decomposition products:** Not determined.

Article 11 - Toxicological Information

- **Acute toxicity:** Not available.
- **LD/LC50:** Not available.
- **Skin corrosion/irritation:** Not available.
- **Serious eye damage/eye irritation:** Not available.
- **Respiratory or skin sensitization:** Not available.
- **Germ cell mutagenicity:** Not available.
- **Carcinogenicity:** No components are listed in IARC, or NTP, or OSHA, or ACGIH.
- **Reproductive toxicity:** Not available.
- **Teratogenicity:** Not available.
- **Specific target organ toxicity - single exposure/ - repeated exposure (GHS):** Not available.
- **Aspiration hazard:** Not available.
- **Potential health effects:**
Inhalation: No data available
Ingestion: No data available
Skin: No data available
Eyes: No data available
- **Signs and Symptoms of Exposure:** No data available
- **Synergistic effects:** Not available.

Article 12 - Ecological Information

- **Eco-toxicity:** No data available.
- **Biodegradability:** Not applicable.
- **Bio-accumulative potential:** Not applicable.
- **Mobility in soil:** Not applicable.
- **PBT and vPvB assessment:** Not applicable.
- **Other adverse effects:** Not applicable.

To place your order, please contact us by phone 1-(604)-232-4600, fax 1-604-232-4601 or by email: orders@signalchem.com or IVD@signalchem.com - www.signalchem.com

FOR IN VITRO RESEARCH PURPOSES ONLY. NOT INTENDED FOR USE IN HUMAN OR ANIMALS.

SAFETY DATA SHEET

Article 13 - Disposal Considerations

- **Disposal methods:** In accordance to applicable national, regional, or local laws and regulations. For additional handling information and protection of employees please refer to Article 7 and 8.
- **Contaminated packaging:** Disposal should be made in accordance to official regulations. Use water or cleansing agents to clean the area.

Article 14 - Transport Information

- **DOT:** Not dangerous goods.
- **IMDG:** Not dangerous goods.
- **IATA:** Not dangerous goods.

Article 15 - Regulatory Information

- **WHMIS Classification:** Non-hazardous.
- **GHS label elements:** Not applicable.
- **Signal word:** Not applicable.
- **Hazard statements:** Not applicable.

Article 16 - Other Information

The above information is believed to be correct but does not purport to be all-inclusive and shall be used only as a guide. SignalChem shall not be held liable for any damage resulting from handling or from contact with the above product. See the Technical Specification, Packing Slip, Invoice, and Product Catalog for additional terms and conditions of sale.

To place your order, please contact us by phone 1-(604)-232-4600, fax 1-604-232-4601
or by email: orders@signalchem.com or IVD@signalchem.com - www.signalchem.com

FOR IN VITRO RESEARCH PURPOSES ONLY. NOT INTENDED FOR USE IN HUMAN OR ANIMALS.