

INSTRUCTIONS FOR USE

Product Name: RNase R, GMP-Grade (20 U/ μ L)**Catalog # RNR01-E59C**

Step 1. In a sterile microfuge tube, add the following components:

2 μ L	10x RNase R Reaction Buffer
1 μ g	RNA Sample
2-4 U	RNase R (20 U/ μ l)
Up to 20 μ L	DEPC H ₂ O

Step 2. Incubate the reaction at 37°C for 10 - 30 min.

Step 3. To inactivate the reaction, incubate at 70°C for 10 min.